

# TEST PROCEDURE



**1**

Fill the test tube (0106) to the 10 mL line with sample water.



**2**

Add one \*Phosphate LR Testab (6836A).



**3**

Use the Tablet Crusher (0175) to crush the tablet.



**4**

Cap (28912) the tube & mix until the tablet has disintegrated.




**5**

Wait 6 minutes. Invert the tube once to mix. Remove cap.



**6**

Place the bottom of the tube inside the dotted lines on the color chart.

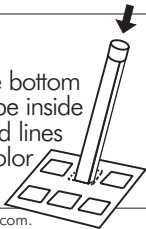

Look down through the tube to match the sample color to a color standard. Record the result as ppb Phosphate.

Store the kit at room temperature.

**LaMotte Company**

PO Box 329

Chestertown, MD 21620

800-344-3100

410-778-3100 (Outside USA)

[www.lamotte.com](http://www.lamotte.com)

\*Potential health hazard. Read MSDS at [www.lamotte.com](http://www.lamotte.com).



low range  
PHOSPHATE  
TEST KIT

Code 3466